Moku:Pro Application Note September 1, 2021

## Laser Frequency Offset Locking

# An FPGA-based, 4-channel Phasemeter for optical phase-locked loop (OPLL)

Stability is vital for sensitive measurement systems. It is the key parameter that determines the precision and accuracy of such systems. Like a voltage reference in a voltmeter, the frequency and phase of a laser must be referenced to a stable source. In this technical note we look at offset phase locking in optical systems as a method to transfer the stability of one optical system to another.



#### Introduction

Optical phase-locking is a common technique for transferring the frequency and phase characteristics of one laser beam to another. It is commonly used in heterodyne metrology, free-space optical communications and spectroscopic applications. In this technical note we discuss the implementation of offset phase locking using a digital phasemeter and characterize the stability of phase locking two lasers.

#### **Optical Offset Phase Locking (in a nutshell)**

To put it simply, offset phase locking stabilizes phase difference between two lasers. This is done by first measuring the phase difference of the lasers and second, by feeding back to modify the frequency of one of the lasers to stabilize the phase difference.

Measuring the phase difference between two lasers is a relatively simple process combining the outputs of the two lasers (via a beamsplitter or similar) and illuminating a photodetector with the combined beam. The result is analogous to a mixing process, and produces an oscillating signal at the difference frequency of the two lasers. We can refer to this as a beat note.

The power at the photodetector can be described through the following equations:

$$P_{PD} = |E_{PD}|^2 = |E_1 + E_2|^2 \tag{1}$$

 $P_{PD}$  and  $E_{PD}$  are the power and electric field at the detector.  $E_1$  and  $E_2$  are the fields of the output of the respective lasers and are given by:

$$E_1 = A_1 e^{-i(\omega_1 t + \phi_1)} \tag{2}$$

$$E_2 = A_2 e^{-i(\omega_2 t + \phi_2)} \tag{3}$$

Where  $\omega_1$  and  $\omega_2$  denote the frequency, and  $\phi_1$  and  $\phi_2$  denote the phase of each laser. Substituting eq (2) and eq (3) into eq (1) we get

$$P_{PD} = \cos((\omega_1 - \omega_2)t + \phi_1 - \phi_2) + higher order terms$$
 (4)

Note that the high order terms are generally outside the bandwidth of the photodetectors. It is important to recognize that even though the beat note contains phase information of the lasers, this information is contained in the argument of the signal and is relatively difficult to use in feedback systems in this form. To extract the phase from the beat note we use a phase detector. A simple phase detector consists of a mixer followed by a low pass filter to transform the information to baseband so that it can be used in a feedback system.

Once at baseband the phase signal can be fed back to one of the lasers to remove any differences between the two lasers. The complete set up is summarized below in Figure 1.


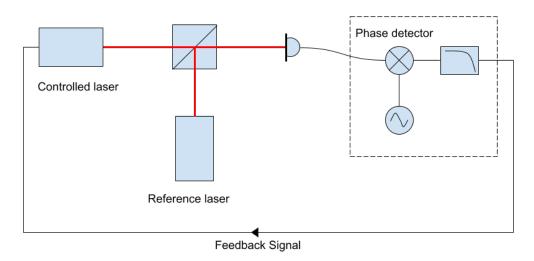


Figure 1: Schematic of a typical offset laser frequency lock system.

#### Phase-locked Loop (PLL) - a different type of phase detector

Although a mixer and low pass filter are adequate for most offset phase locking system, they are not without their limitations. For one, the range of the mixer-filter combination is limited to  $\pm \pi/2$ , and the phase output from the system is linear only when very close to zero. These range and linearity issues often make it difficult to deal with systems with large fluctuations. Using an additional phase-locked loop (PLL), instead of a standard mixer filter, for phase detection may help in these situations.

PLL is a technique used to establish a stable frequency and phase relation between two oscillators. It is widely used in modern electronics, integrated chips, and many other fields. A PLL has three main components: a phase detector, a loop filter, and a controllable/tunable oscillator. The phase detector generates a signal that contains a DC component that is (near) proportional to the phase difference between two signals. This signal is then sent to a loop filter (lowpass and/or PID) to remove the higher harmonics from the phase detector. The filtered error signal is then sent to an oscillator where the output frequency is controlled by the input DC voltage. By forming this closed loop and tuning the loop filter, a stable phase relation of the two oscillators can be achieved.

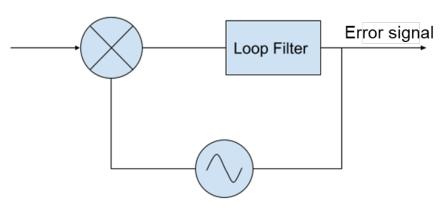

Tunable oscillator

Figure 2: A block diagram of a typical PLL.

The PLL can assist in initial lock acquisition, removal of non-linear affects (such as cycle slipping), and ensure a more robust lock.

## Offset Phase Locking with Moku:Pro

We demonstrated Moku:Pro's Phasemeter offset phase locking architecture using two nonplanar ring oscillators (NPRO) lasers. The beams of the master and slave lasers were combined by a beamsplitter and interfered at a photodiode as shown in Figure 3. The beat note was connected to Moku:Pro's input 1. The feedback signal was then connected to the frequency controller of the following laser.

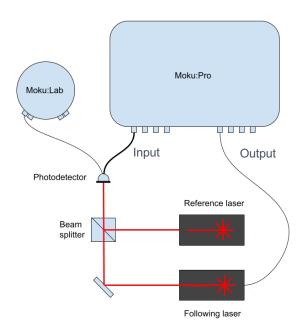


Figure 3: Instrument setup for laser frequency offset locking.

### Setting up the Phase Lock

Before the phase lock can be implemented the system needs to be tuned to operating range. To acquire lock, the thermal actuators were used to coarsely tune the lasers to produce a beat note within the 600 MHz bandwidth of the Moku:Pro inputs. Once in range, the Phasemeter can be set to track the beat note by either using the auto-acquire function, or by manually setting the frequency. Detailed information about the Phasemeter can be found in Ref [5], and the Phasemeter User Manual.

Set the outputs to use phase and select a voltage scaling (this can be treated as gain in a typical control loop). You can generally start off with a small gain a gradually increase the gain to optimize the system.

Moku:Pro Application Note September 1, 2021



Figure 4: Control loop initial gain can be set under the 'Output' pane via 'Scaling'.

In manual operation the initial frequency can be used to tune the frequency of the offset phase lock to the required frequency.

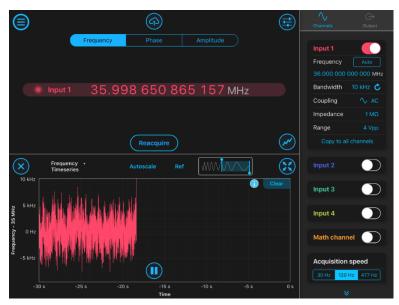


Figure 5: The offset frequency can be tuned via the "Frequency" under "Channel" pane.

#### Lock Performance

The performance of the phase lock was measured using a separate Phasemeter running on a <u>Moku:Lab</u>. Figure 6 shows the locked and free running frequency (a), and phase (b) over a measurement of 60 seconds. Both phase and frequency fluctuations between the two lasers can be clearly seen to be significantly reduced.

Moku:Pro Application Note September 1, 2021

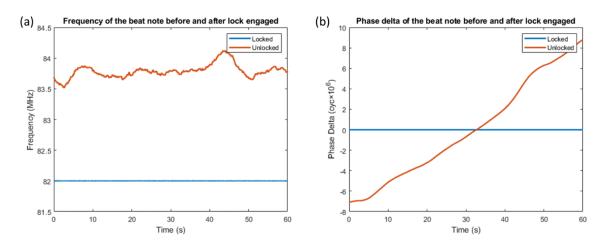


Figure 6: Frequency (a) and phase (b) of the offset beat note locked and unlocked.

Taking the amplitude spectral density of the frequency, we see that an improvement in the stability of over 4 orders of magnitude was measured, with relative frequency stability of 1 Hz/ $\sqrt{\text{Hz}}$  measured down to 0.1 Hz.

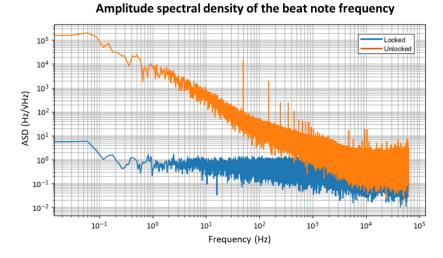


Figure 7: Amplitude spectral density of the offset beat note before and after the lock was engaged.

#### Conclusion

Laser offset locking maintains the frequency difference between a master and a slave laser. The dynamic range for the phase error signal in such a system is typically beyond  $2\pi$ , where a mixer type phase detector may fail to continuously track the signal. Moku:Pro's Phasemeter implements four independent PLL phase detectors with automatic phase unwrap functions. A stable lock between two lasers with an 82 MHz offset was demonstrated with a frequency stability better than  $1 \text{ Hz}/\sqrt{\text{Hz}}$  at 10 Hz.

### Acknowledgment

We would like to thank The Australian National University for providing us the details of their experiment, explanation of use of Moku:Pro and feedback. The experiment at the ANU was supported by the ARC Centre of Excellence for Gravitational Wave Discovery.

#### References

[1] Schünemann, U., H. Engler, R. Grimm, M. Weidemüller, and M. Zielonkowski. "Simple scheme for tunable frequency offset locking of two lasers." *Review of Scientific Instruments* 70, no. 1 (1999): 242-243.

- [2] Bell, S. C., D. M. Heywood, J. D. White, J. D. Close, and R. E. Scholten. "Laser frequency offset locking using electromagnetically induced transparency." Applied physics letters 90, no. 17 (2007): 171120.
- [3] Thorpe, James I., K. Numata, and J. Livas. "Laser frequency stabilization and control through offset sideband locking to optical cavities." *Optics express* 16, no. 20 (2008): 15980-15990.
- [4] Hsieh, Guan-Chyun, and James C. Hung. "Phase-locked loop techniques. A survey." *IEEE Transactions on industrial electronics* 43, no. 6 (1996): 609-615.
- [5] Shaddock, D., Ware, B., Halverson, P. G., Spero, R. E., & Klipstein, B. (2006, November). Overview of the LISA Phasemeter. In AIP conference proceedings (Vol. 873, No. 1, pp. 654-660). *American Institute of Physics*

### Questions or comments?

Please contact us at <a href="mailto:support@liquidinstruments.com">support@liquidinstruments.com</a>